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# Automatic film applicator coater

## ZAA 2300

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Manual applications may have irregularities caused by uneven pressure or non-uniform drawing speed. The ZAA 2300 Automatic film applicator coater ensures an accurate and reproducible coating.

## Features

- double sided glass plate for dual-purpose operation: one side with printing blanket for wire-bar applicators and profile rods, on the other side glass surface for other applications
- option: vacuumplate for substrates; the vacuum created fixes the substrate in place evenly

## Standard delivery

- 1 automatic film applicator coater
- 1 reversible, double sided glass plate / printing blanket
- 1 spirit level
- 1 certificate of manufacturer

## Options

- ACC121 precision vacuumplate
- ACC039/ACC122 (230V/115V) vacuum pump with hose
- ACC225.230/ACC225.115 (230V/115V) heating plate / ACC188/ACC189 (230V/115V) heated precision vacuum plate all for temperature up to 150°C (302°F), those options can only be used together with a isolation plate equipped automatic film applicator coater (2300H/FH/FFH)
- ZTC 2200 Test charts
- film applicators with maximum outer width of 300 mm (11.81")

## Handling

- choose the applicator(s)
- choose the appropriate base (glass surface, printing blanket for wire bar applicators and profile rods or vacuumplate)
- mount wire bar applicator or profile rod to the ZAA 2300 Automatic film applicator coater and fix with the weight support
- position other applicators in front of the weight support
- adjust the desired application length
- fix the desired substrate, e.g. test chart, with the start position clamp
- set the desired drawing speed
- pour the product to be tested in front of the applicator in drawing direction
- start the ZAA 2300 Automatic film applicator coater
- remove the applicator, the weight and the substrate
- afterwards clean the applicator and, if necessary, the base